



KERN & Sohn GmbH

Ziegelei 1

D-72336 Balingen

E-mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: www.kern-sohn.com

Printer configuration

KERN YKH-01

Version 1.0

2025-03

GB



TYKH-01-B-ZB_KFG-e-2510

KERN YKH-01

Please refer to the operating instructions supplied with the printer for the relevant printer settings.



KERN YKH-01

Printer settings and scale settings must match for proper data transfer.
The enclosed serial adapter must also be connected.

Standard printer settings

Interface	RS232
Baud rate	38400
Data bit	8 bit
Stop bit	1 bit
Parity	None
Handshaking	DTR/DSR

Type Scale series	Interface cable	Comment
TPFB-B; TEWJ-B	CFS-A01	
EMB-V; PCD; 440; EN-D/A	572-926	
TPCB-A/TPCJ-A; TFCB-A/-B; TFKB-B, TCKE-B; TCKE-A (small); TIFC-C; TBFC-C; TUFC-C; TNFC-C; TKFC-TM-C; TDS-A; TCDS-A; TPKS-A/TPKJ-A; T572-A	KUP-01	
TIXC-A; TSXC-A; TBXC-A; TUXC-A; TNXC-A; TKXC-TM-A	KUM-01	Install interface: KUM-01
TALSG-A/TALJG-A; TPLSG-ATPLJG-A	572-926	
PNS; PNJ	MLB-A05	
EW-N; EG-AM	474-926	
TPWS-A	MLB-A05/PWS-A02	
PBS; PBJ-N	770-926	
TPES-B/TPEJ-B	474-926/MLB-A05	
TFES-A/TFEJ-A	MLB-A05	
ADB-A; TADB-B; TADJ-A/-B	MLB-A05	
TAXS-A/TAXE-A	MLB-A05	
TADS-A/TADT-A; TPDS-A/TPDT-A	572-926	
TACS-A/TACJ-A	MLB-A05	
ABT-NM	770-926	
TABP-A	MLB-A05	
DAB/TDAB-A	MLB-A05	
DBS	MLB-A05	
TDLG-A	572-926	

Type Scale series	Interface cable	Comment/ DIRECT
MPS-NM/MWS-NM/MCB-NM	MPS-A08	
TMPN-A / TMWN-A / TMCN-A	KUP-01	
MPE/MTA-NM/TMPE-EM-A	CFS-A01	
MPC-NM/MPC-LM	CFS-A01	Install interface: MPC-A01
MWA-NM	CFS-A01	
CPB-N/CPB-DM	CFS-A01	
CCA	CCA-A01	
CCS	CFS-A01	
TFCD-A; TCPE-A; TEFC-A	CFS-A01	
RFC, TRFC-A	CFS-A01	
TSFB-A; KFN-TM	CFS-A01	Install interface: KFN-A01
IOC-M, TIOC-A, BID-M, BID-D, TBID-A, UID-M, UID-DM, KIB-TM	EOC-A12	